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Baoding Longer Precision Pump Co., Ltd.

⚠ Important Information:

Please read operating manual carefully before operation.

⚠ Warning:

- Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment. Please check the tubing frequently and change the tubing in time.


Instruction

YZII series pump heads have good appearance and compact size. The tubing retention set is a linkage, it is convenient to load the tubing; Accept several tubing sizes for a wide flow range; The rollers adopt high Performance materials and has good wearing property; The housing material is PESU. It has perfect rigidity and structure property. It is stable and autoclavable.

Type	Number of Rollers	Roller Material	Housing Material	Weight (kg)
YZ II 15	3	RULON [®] J	PESU	0.35
YZ II 25				

Type	Applicable Tubing				Flow Rates (mL/min)
	Tubing Type	ID×Wall Thickness(mm)	OD (mm)	Speed (rpm)	
YZ II 15	13 [#]	0.8 × 1.6	4.0	1-600	0.06-42
	14 [#]	1.6 × 1.6	4.8		0.27-162
	19 [#]	2.4 × 1.6	5.6		0.41-250
	16 [#]	3.2 × 1.6	6.4		0.77-466
	25 [#]	4.8 × 1.6	8.0		1.6-980
	17 [#]	6.4 × 1.6	9.6		2.5-1500
	18 [#]	8.0 × 1.6	11.2		3.8-2280
YZ II 25	15 [#]	4.8 × 2.5	9.8		1.7-966
	24 [#]	6.4 × 2.5	11.4		2.8-1653
	35 [#]	7.9 × 2.5	12.9		4.1-2495
	36 [#]	9.6 × 2.5	14.6		5.3-3000

Tubing Loading

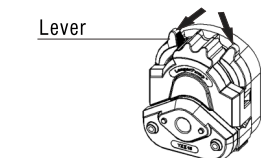
1. Pull the **Levers** according to the arrow head directions simultaneously and open the **Compression Block**.


Prompt


The levers in both sides need to be pulled simultaneously.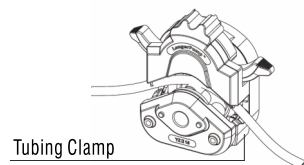

2. The **Tubing Clamp** and the **Compression Block** is linkage. Tubing retention is automatic.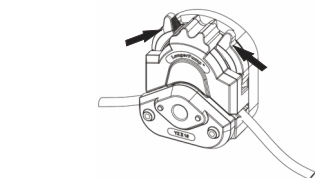

Structure

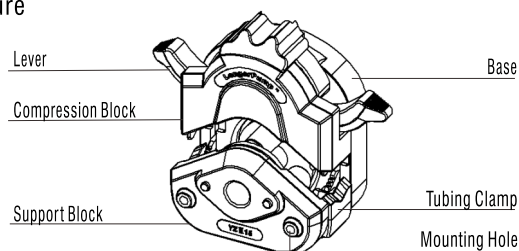The material of pump head is not organic solvent resistant.

Maintenance

- When the pump is idle, we recommend you to release the tubing from pressure. This helps to protect the tubing from unnecessary strain and prolongs its service life
- Keep rollers clean and dry. This will prolong the service lives of tubing and pump head.
- The materials of pump head are not organic solvent and aggressive liquids resistant. Please pay attention when using.


If a trouble happens, please contact us or our dealers.


Warranty

The warranty period for this product is one year. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

Pump Head Installation

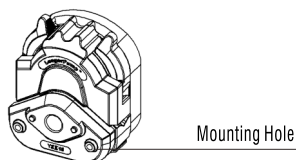
- Insert the two mounting screws into the mounting holes of the pump head. Insert the flat end of the pump head's shaft to the slot of drive's coupling. Then tighten the mounting screws to connecting hole of drive. Stack pump heads and long mounting screws.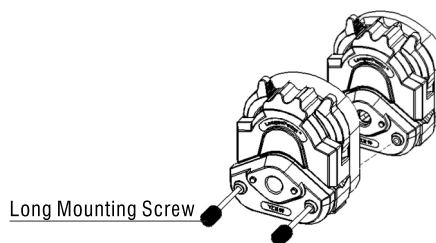

Try your best to tighten the two screws in the same degree. Do not over-tighten the screws to prevent the support block from deforming and noises during running.

Dimensions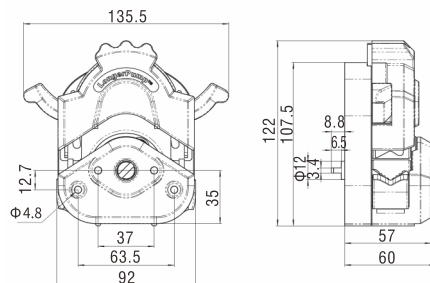

Unit: mm